User Guide

Elcometer 2350, 2351,
2352, 2353 & 2354

Viscosity Flow Cups

CONTENTS

en	1	Overview
	2	Box Contents
	3	Test Procedure
	4	Care & Maintenance
	5	Spares & Accessories
	6	Warranty Statement
	7	Technical Specification

For the avoidance of doubt, please refer to the original English language version.

Elcometer Viscosity Flow Cups are packed in a cardboard package. Please ensure that this packaging is disposed of in an environmentally sensitive manner. Consult your local Environmental Authority for further guidance.

elcometer® is a registered trademark of Elcometer Limited, Edge Lane, Manchester, M43 6BU.
United Kingdom

All other trademarks acknowledged.

© Elcometer Limited 2009-2021. All rights reserved. No part of this document may be reproduced, transmitted, transcribed, stored (in a retrieval system or otherwise) or translated into any language, in any form or by any means (electronic, mechanical, magnetic, optical, manual or otherwise) without the prior written permission of Elcometer Limited.

1 OVERVIEW

Elcometer Viscosity Flow Cups are easy-to-use instruments for measuring the consistency of liquids.

The cup is supported in a stand^a and is filled with the liquid under test. The time taken for the liquid to drain through the orifice in the bottom of the cup is measured.

The measured kinematic viscosity is generally expressed in seconds (s) flow time, which can be converted to Centistokes (cSt), if stipulated in the Standard. The Elcometer 2400 Viscosity Conversion Disc^a, Elcometer ElcoCalc™ Mobile App (see www.elcometer.com) or the look-up tables in Appendix 1 (after page jp-7) can be used for this purpose.

A wide range of cups with different orifices is available for measurements between 5 cSt and 5100 cSt.

2 BOX CONTENTS

- Elcometer Viscosity Flow Cup
- Batch or Cup Specific Calibration Certificate^b
- User Guide

3 TEST PROCEDURE

3.1 BEFORE YOU START

- 1 Select a cup which gives a flow time of between 30 seconds and 100 seconds.
- 2 Ensure the cup and the orifice are clean and free of debris.
- 3 The liquid being tested must be homogeneous and must not contain any bubbles. Use freshly strained samples only.
- 4 The liquid being tested must be newtonian. To determine whether a liquid is newtonian, carry out the following measurements using the instructions given in Section 3.2 'Test Procedure' on page en-3:
 - Fill the cup and measure flow time immediately after filling.
 - Fill the cup, wait one minute, then measure flow time.

If the difference between the two flow times is greater than 10%, the liquid is considered to be non-newtonian and cannot be tested using the Viscosity Cup.

- 5 Measure and record the temperature of the liquid.

^a Available to purchase as an optional accessory, see Section 5 'Spares & Accessories' on page en-5.

^b Depending on part number ordered, see Section 7 'Technical Specification' on page en-6 for details.

3 TEST PROCEDURE (continued)

en

3.2 TEST PROCEDURE

- 1 Place the cup into the stand^a and level the cup using the glass plate and bubble level supplied with the stand (Figure 1).
- 2 Allow the temperature of the cup and the sample to stabilise at the agreed temperature. Check the temperature of the sample before filling^c.
- 3 Place a finger over the cup orifice to close the orifice (Figure 2).
- 4 Pour the liquid gently into the cup, avoiding the formation of air bubbles (Figure 2).
- 5 Slide the glass plate over the rim of the cup to remove excess sample. Avoid the formation of air bubbles between the glass plate and the liquid (Figure 3).
- 6 Remove your finger from cup orifice and wait for a few minutes to allow any air bubbles to rise to the top of the liquid.
- 7 Remove the glass plate and start the stopwatch^a (Figure 4).
- 8 Watch the flow of liquid from the orifice and stop the stopwatch^a the moment the flow breaks for the very first time (Figure 5).
- 9 Repeat the test.
- 10 If the test results do not differ by more than 5%, calculate the average of the two test results and record the result.

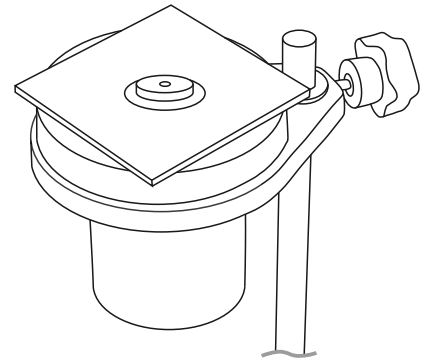
Figure 1

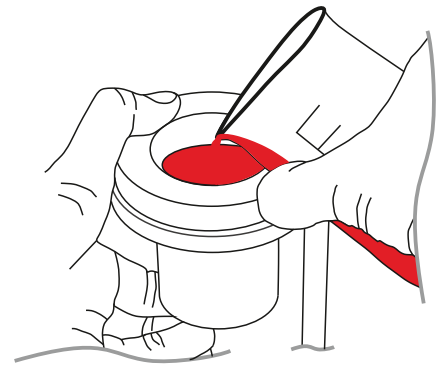
Figure 2

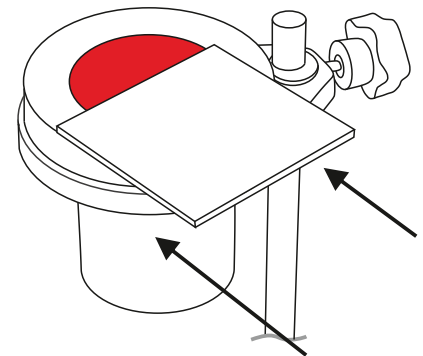
Figure 3

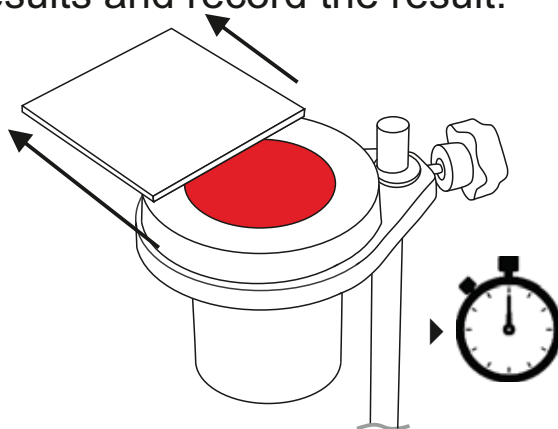
Figure 4

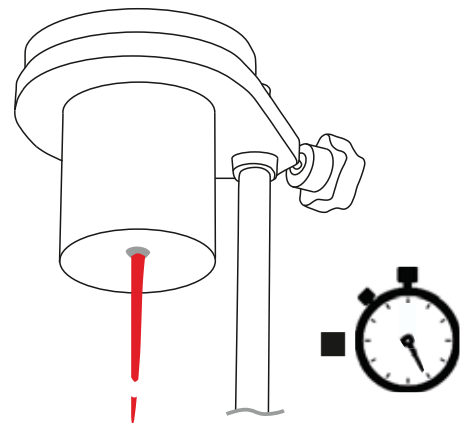
Figure 5

^a Available to purchase as an optional accessory, see Section 5 'Spares & Accessories' on page en-5.

^c If the test environment is not temperature stabilised a double-walled thermo-jacket can be used, see Section 5 'Spares & Accessories' on page en-5.

3 TEST PROCEDURE (continued)

The measured kinematic viscosity is generally expressed in seconds (s) flow time, which can be converted to Centistokes (cSt), if stipulated in the Standard. The Elcometer 2400 Viscosity Conversion Disc^a, Elcometer ElcoCalc™ Mobile App (see www.elcometer.com) or the look-up tables in Appendix 1 (after page jp-7) can be used for this purpose.

3.3 AFTER TEST

Clean the viscosity cup and all the equipment used using a suitable solvent.

DO NOT use wire brushes, metal scrapers, metal files or other metallic tools for cleaning.

After cleaning, ensure that all materials are removed and that the instrument is dry.

4 CARE & MAINTENANCE

Elcometer Viscosity Flow Cups are designed to give many years reliable service under normal operating conditions.

- Regularly inspect the cup for damage to the orifice, scratches on the internal surface or deformation as such flaws will affect the readings and the cup may have to be replaced.
- To check for wear, use Elcometer 2410 Viscosity Cup Standard Calibration Oils^a in place of your liquid and measure the flow time. If wear is detected, contact Elcometer or your local Elcometer supplier.
- Regular calibration checks over the life of the cup are a requirement of quality management procedures e.g. ISO 9000 and other standards.

^a Available to purchase as an optional accessory, see Section 5 'Spares & Accessories' on page en-5.

5 SPARES & ACCESSORIES

en 5.1 VISCOSITY CUP STANDS

Using a viscosity cup stand ensures that the cup is positioned correctly to carry out the test.

Each stand is supplied with bubble level; to ensure the cup is parallel to the surface and glass draw down plate; to retain the test sample until the user is ready to begin the test. The glass draw down plate also provides a surface for the bubble level.

A stand complete with thermo jacket is also available for heating test samples if the temperature of the test environment is unstable.

Description

Viscosity Cup Stand

Viscosity Cup Precision Stand

Viscosity Cup Precision Stand with Thermo jacket

Bubble Level

Glass Draw Down Plate

Thermo jacket

Part Number

KT002400N201

KT002400N001

KT002400N202

KT002400P001

KT002400P999

KT002400N002

5.2 STOPWATCH

To time the flow of the liquid from the viscosity cup.

Description

Elcometer 7300 High Precision Stopwatch

Part Number

K0007300M201

5.3 VISCOSITY CONVERSION DISC

The measured kinematic viscosity is generally expressed in seconds (s) flow time. This can be converted to Centistokes (cSt) using the Elcometer 2400 Viscosity Conversion Disc.

Description

Elcometer 2400 Conversion Disc

Front: No.4 cups according to AFNOR, BS, NF, ASTM, DIN, Zahn 2

Back: No.3-4-5-6 cups according to ISO and Zahn 3

Part Number

KT002400N003

5 SPARES & ACCESSORIES (continued)

5.4 ELCOMETER 2410 VISCOSITY CUP STANDARD CALIBRATION OILS

In order to check the viscosity cup's calibration or to certify it for ISO purposes, it is imperative that viscosity cup standard calibration oils are used.

Supplied in ½ litre (1 pint) bottles complete with calibration certificate, standard oils have a specific drain time, dependent upon the viscosity cup type (Ford, DIN, ISO etc.) and the orifice size or cup number used.

Specific calibration oils can only be used with specific flow cups - refer to the table below to determine which calibration oil is required for each cup.

Viscosity Oil Part Number	Flow Cup Type / Number			Kinematic Viscosity at 25°C (77°F) ^e
	DIN	ASTM / FORD	ISO	
K0002410M021	-	2	3	34cSt
K0002410M022	4	3	4	120cSt
K0002410M023	4	4	6	230cSt
K0002410M024	4	-	6	460cSt

Material Safety Data Sheets for the Elcometer 2410 Viscosity Cup Standard Calibration Oils can be downloaded via our website:

http://www.elcometer.com/images/stories/MSDS/elcometer_2410m021.pdf

http://www.elcometer.com/images/stories/MSDS/elcometer_2410m022_m023.pdf

http://www.elcometer.com/images/stories/MSDS/elcometer_2410m024_m025_m026.pdf

6 WARRANTY STATEMENT

Elcometer Viscosity Flow Cups are supplied with a one year warranty against manufacturing defects, excluding contamination and wear.

7 TECHNICAL SPECIFICATION

Body Material	Anodised Aluminium	
Orifice Material	Stainless Steel	
Accuracy	Elcometer 2350:	3%
	Elcometer 2351:	10%
	Elcometer 2353:	3%
	Elcometer 2354:	2.5%

^e Nominal values.

7 TECHNICAL SPECIFICATION (continued)

en

Part Number [‡]	Cup Type / Number	Orifice Diameter (mm)	Applicable Standards	Range (cSt)	Certificates Available [‡]	
					Batch	Cup
K0002350M001	DIN 2	2.00	DIN 53211 (Cup 4 only)	-		○
K0002350M002	DIN 4	4.00		96 - 683	◆	□
K0002350M003	DIN 6	6.00		-		○
K0002350M004	DIN 8	8.00		-		○
K0002351M001	FORD / ASTM 1	1.90	ASTM D1200 (All cups)	10 - 35	◆	□
K0002351M002	FORD / ASTM 2	2.53		25 - 120	◆	□
K0002351M003	FORD / ASTM 3	3.40		49 - 220	◆	□
K0002351M004	FORD / ASTM 4	4.12		70 - 370	◆	□
K0002351M005	FORD / ASTM 5	5.20		200 - 1200	◆	□
K0002352M001	AFNOR 2.5	2.46	NF T30-014 (All Cups)	-		○
K0002352M002	AFNOR 4	4.00		-		○
K0002352M003	AFNOR 6	6.00		-		○
K0002353M001	ISO 3	3.00	ISO 2431 [†] (Cups 3, 4, 5 & 6) ASTM D 5125 (Cups 3, 4, 5 & 6)	7 - 42	◆	□
K0002353M002	ISO 4	4.00		34 - 135	◆	□
K0002353M003	ISO 5	5.00		91 - 326	◆	□
K0002353M004	ISO 6	6.00		188 - 684	◆	□
K0002353M005	ISO 8	8.00		-		○
K0002354M003	BS 4	3.97	BS 3900 A6:1971 (Cup 4 & 5)	89 - 340	◆	□
K0002354M004	BS 5	4.76	AS/NZS 1580.214.2 (Cup 4 only)	79 - 441	◆	□

- ◆ Batch Certificate supplied as standard.
- Optional Calibration Certificate Available: Efflux Time Certificate
- Optional Calibration Certificate Available: Dimensional Certificate

[‡] Add 'C' to the end of the sales part number for cup supplied complete with Efflux Time or Dimensional Calibration Certificate as indicated. Calibration Certificates must be requested at time of order, they can not be supplied retrospectively.

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© Elcometer Limited 2009-2021. All rights reserved. この文書の一部または全部を、Elcometer Limitedの事前の書面による許可なく、いかなる形式や方法（電子的、機械的、磁氣的、工学的、手動を問わず）によっても、複製、転送、保管（検索可能なシステムかどうかを問わず）、または他の言語に翻訳することを禁じます。